


# FullRanger 100bp DNA Ladder

(100bp - 5,000bp)

Catalogue Number: 11800

100 gel lanes

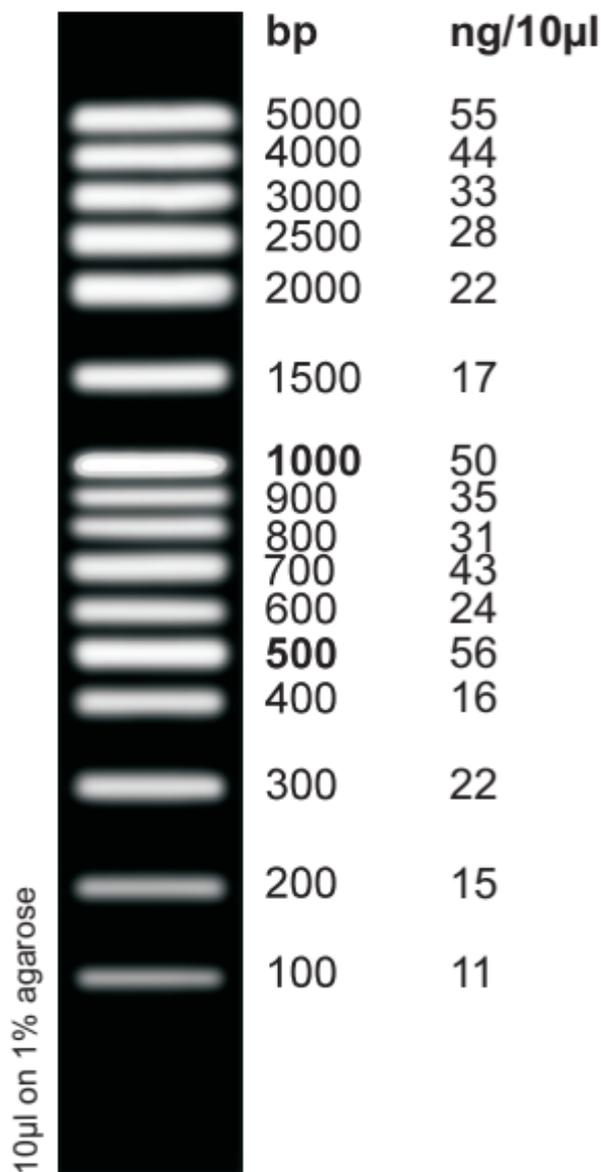



## **FullRanger 100bp DNA Ladder (100bp - 5,000bp)**

Catalogue Number: 11800 (100 gel lanes)

### **Description:**

High quality ready-to-use double-stranded DNA ladder derived by recombinant technology for precise on-gel size and mass determination.

### **Features:**

- Precise band sizes and masses for densitometric quantification.
- Sixteen discrete bands, ranging from 100bp to 5,000bp.
- Higher intensity band at 500bp and 1,000bp for easy reference.

### **Contents:**

1mL of premixed DNA ladder (0.5µg/10µL) in loading buffer (10mM EDTA, 10% glycerol, 0.015% bromophenol blue, and 0.17% SDS).

### **Instructions:**

Mix thoroughly. For best results, load 10µL of DNA ladder per well. For precise mass determination with a densitometer, stain gel after electrophoresis using 0.5µg/mL ethidium bromide for 30-40 minutes. The accompanying diagram shows the size and mass for each band based on 10µL ladder per well.

### **Storage:**

Stable at room temperature. For longer term storage, -20°C is recommended.

This ladder was standardized using 10µL of DNA per lane on a 0.8 cm thick, 13 x 15 cm, 1.0% agarose gel run in TAE buffer.

Tech support: [techsupport@norgenbiotek.com](mailto:techsupport@norgenbiotek.com)

Orders: [orders@norgenbiotek.com](mailto:orders@norgenbiotek.com)

Toll Free: 1-866-667-4362

Fax: 905-227-1061

[www.norgenbiotek.com](http://www.norgenbiotek.com)