
cf-DNA/cf-RNA Preservative Tubes

Mixing Instructions

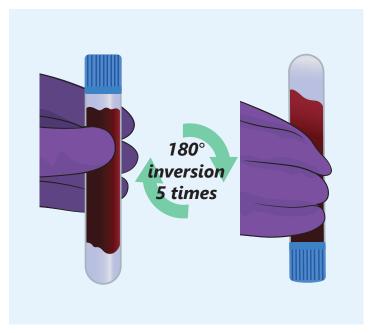



Norgen Biotek Corp. 3430 Schmon Pkwy, Thorold, ON, Canada, L2V 4Y6 www.norgenbiotek.com | info@norgenbiotek.com

IMPORTANT!

Immediately after blood is drawn into the tube, invert the tube five times (x5) to ensure that the preservative makes uniform contact with the sample.

One inversion is a complete turn of the wrist 180 degrees and back.

DO NOT MIX the contents of the cf-DNA/ cf-RNA Preservative Tube again once this step is complete.

Cat. 63950, 63960